



RSolv[®]
CHROMATOGRAPHY PRODUCTS

LAB CLEAN

Best Solution For Your Laboratory Glassware



Lab Clean generally used in Research centers,
Path labs, Diagnostic centers, Hospital and many more.

Ingredients :

LABSA/DBSA -2%, SODA ASH -2%, STP -20%,
SLES -5% Aqua Solution -81%

Method of Use :

Pour 70 ml of Lab Clean in a bucket and
drop all glassware/plasticware. Take out after 15min
and wash with regular water



Mfg. Date	: Aug 2023
Expiry Date	: One year from the date of mfg
Batch No.	: R-008
ML No.	: RAK-2721
M.R.P.	: 1500.00

Distributed By:

RIKI GLOBAL TRADING PVT LTD

Ph.No : 040 - 42217787, 9908246695

Email Id : info@rikiglobal.com

Website : www.rikiglobal.com